

Heating Mantle

USER MANUAL

I. Introduction

The instrument is designed for heating liquids in schools, laboratories or factories. This device is not suitable for using in residential areas.

II. Operating

- Place the equipment on level worktable, and then put the container with liquid inside into heating mantle.
- Switch on the power accords with the machine, then power indicator light will be lighten; turn on the power of regulation knob, and turn the knob clockwise, then the working indicator light will be lighten; in the process of regulating, the light intensity changes according to different regulation position, and the temperature rises as well.

Notes:



- The power must accord with the machine.
 - Make sure the power line has safety distance from the heating mantle.
 - When the machine meets fault, please cut off the Electricity first.
-

III. Faults

- Instruments can't be power ON
- Check whether the power cable is plugged

If these faults are not resolved, please set the instruments to factory default setting, or take the unit to your technical service center, or contact with the manufacturer.

IV. Maintenance and Cleaning

- Proper maintenance can keep instruments working in a good state and lengthen its lifetime.
- Be careful not spray the cleanser into the instrument when cleaning.
- Unplug the power line when cleaning.

V. Storage and transportation

- Keep it in dry and clean room with good ventilation and no corrosive gas
- prevent it from wetting by the rain and avoid violent collision in transportation.

VI. Working condition

Ambient temperature: 5~40°C;

Ambient humidity: ≤90%;

Voltage: 220V±10%, 50/60Hz